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*FILTRATION AND VACUUM - PUMPS*

*FILTRATION ET VIDE - POMPES*

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### FILTRATION AND VACUUM DESICCATORS

**Material:** Polypropylene - Polycarbonate

Implosion proof desiccators. Will take vacuum of 740 mm Hg with loss not exceeding 20 mm Hg in 24 hours. The transparent polycarbonate cover seals onto a neoprene 'O' ring set into the flange of the polypropylene base. A little grease spread on the 'O' ring will assist in establishing initial vacuum. Within the desiccator is a removable pan for holding a drying agent. Vacuum release tap/non return valve assembly is included; specially grooved stopper in this assembly allows air, when required, to be re-admitted very slowly to avoid disturbing the contents.

**Usage Instructions:** remove the cover in PC - insert the samples in the pan - open the vacuum retention valve (PE needle)  
- connect the tube of the pump to the desiccator spigot in order to create vacuum - close the vacuum retention valve

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### DESICCATOR PLATES

**Material:** Polypropylene

These plates made of polypropylene are extremely useful as a support for crucibles, Petri dishes etc. Use at room temperature. Not suitable for use with hot crucible.

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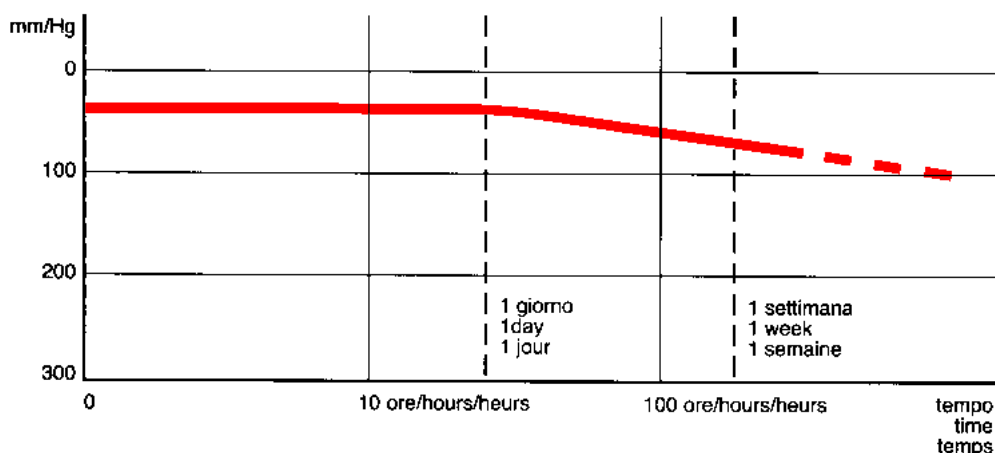
**SPARE VACUUM RETENTION VALVE**

Material: Polycarbonate with Polyethylene Needle

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**WATER JET PUMP**

Material: Polypropylene

This pump will work on pressures up to 10 Kg/cm<sup>2</sup>. Built in non return valve to eliminate back flow. Useful for sucking to waste any split chemicals as well as for filter work. Easily dismantled for cleaning. Autoclavable.

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
The graph below details the performance of the pump under different conditions. Example: Operating pressure of water: 1 Kg/cm<sup>2</sup>. Attained Vacuum: 15 mm/Hg. Will evacuate a 5 litre container in about 20 minutes.

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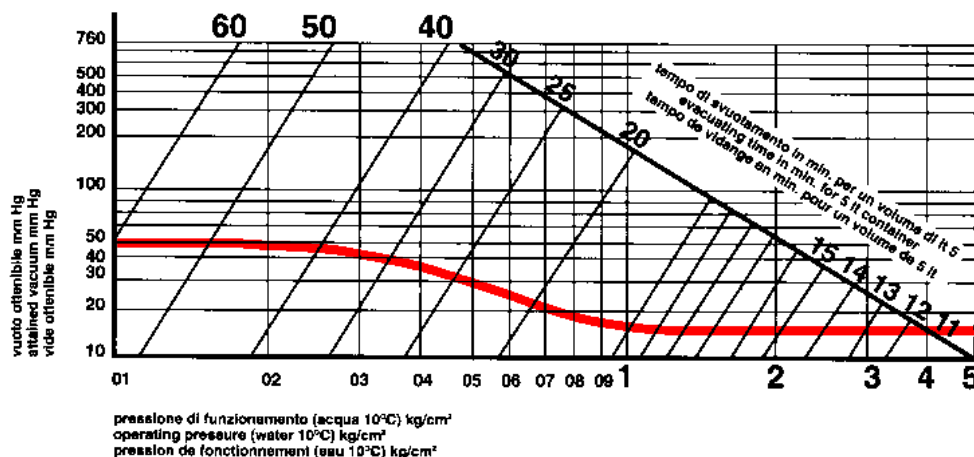**Materiale:** Polypropylene

**Inner Disc Material:** Dutral

Prevents water back-flow when pressure falls. Useful for any tubing system and vacuum line. Maximum safe working pressure: 2 BAR (equal approx. 28.4 P.S.I.). Minimum safe working pressure 0,07 BAR at room temperature (20°C)

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#### FUNNELS, BUCHNER

**Material:** Polypropylene

Robust, lightweight. Unbreakable and resistant to corrosion and heat. Two piece construction. Top and bottom can be separated for easy and total cleaning. Autoclavable. Excellent chemical resistance.

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#### DISCS FOR BUCHNER FUNNELS

**Material:** HDPE

Useful interface between filter paper and perforated Buchner plate. Avoids clogging of filter holes. Texture 0,5 mm.

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#### HAND OPERATED VACUUM PUMP

**Material:** High Impact Polystyrene

Light, portable hand operated vacuum pump. Attains a vacuum of 625 mm Hg (25"). Pumping rate is 15 ml per stroke. Built in vacuum release valve. Can be used as pressure pump by transferring release valve cap to vacuum port. Nozzle fits standard 6 mm (1/4") i.d. tubing.

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**ANALYTICAL FUNNELS**

**Material: Polypropylene**  
 Designed specifically for analytical chemistry. Each size takes a standard size filter paper. Body of funnel is exact 60° angle. Provided with external and internal ribs.  
 Autoclavable. **Suitable for foodstuff.**

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**LONG-STEM FUNNELS**

**Material: Polypropylene**  
 The 60° angles sides are provided with ribs for fast filtration and allow the use of standard size filter papers.  
**Suitable for foodstuff.**

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**SUCTION PUMP FOR STRONG ACIDS**

**Material: PE**  
 Bellows-action siphon pump for high-strength acids. Includes tap.

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**VISCOUS LIQUIDS PUMP**

**Material: Polypropylene**  
 Dispensing pump made of polypropylene operated by a unique lever-action stroke. Delivery of approximately 10 ounces (ab. 300 ml) per stroke, with capability of 6 gallons per minute (ab. 22-23 litres/min). Will handle the most viscous materials (equivalent to SAE 90 oil). Spigot orientation of 360° allows versatility in dispensing. Complete with shank

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### SOLVENT TRANSFER PUMPS

**Material: Heavy nickel chrome plated**

Dispensing pump for solvent transfer with heavy-nickel chrome plated body. Model M45 (Art. N. 2005) fits bottles and buckets with any size and type of openings. Model M855 fits 2" openings in any size drum. NOTE: Solvent pumps are NOT suitable for use with ACID. Shank (Ø2.6 cm) not included with kit.

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